pipetman[®]

M connected

Quick Guide

EN | FR

*Discover the new experience
of connected pipetting on
www.gilsonconnect.com*

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 **GILSON[®]**

Parts Checklist

- PIPETMAN® M Connected pipette
- Power Supply, 5V with AC adapter and cable
- Battery tags (qty 4)
- Ejector tags (qty 4)
- Tip ejector extension (P10M model only)
- PIPETMAN® M Connected Quick Guide
- Safety bag
- Certificate of Conformity with bar-code sticker
- Lubricant (except for the P10M and multichannel models)

Installation

NOTICE

Fully charge the pipette (for approximately three hours) before use.

Use only one of the approved charging devices:

- Power supply, included in box
- Stand adapter, for use with Gilson CARROUSEL™ Stands, sold separately (part number F807023)
- POWER CARROUSEL™ charging pipette stand, sold separately (part number FB1001)

NOTE

The battery can take a few minutes to wake up.

Other References

This Quick Guide provides you with basic information to get started with PIPETMAN M Connected.

Make sure you have read PIPETMAN M Connected User's Guide, available at: www.gilson.com/pipetmanmconnected.

Visit www.gilson.com for additional information.

Get Started

All menu options (refer to [Menu Options](#) on page 10) for PIPETMAN® M Connected can be reached with two buttons.

Push Button

Turn the push button left or right to set the volume. Push to pipette.

Menu Button

Access pipette modes and adjust functions.

Figure 1

Buttons on the PIPETMAN® M Connected

NOTE

After three minutes of inactivity, the pipette display will turn off and enter sleep mode. Press the push button to reactivate the pipette. The last setting and pipetting mode will appear on the screen.

Use the Push Button

- Press the push button to turn on or wake up the startup screen.
- A self-calibration test will begin as the firmware resets the piston.
- The pipette is preset to **PIPET**, also known as standard pipetting.

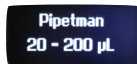
Figure 2
Volume range displayed on startup screen

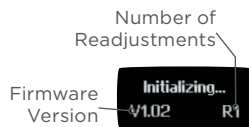
Figure 3
Screen displayed during calibration

Figure 4
Volume screen with pipetting mode displayed

Adjust the Volume

- Quickly turn the push button halfway. 'Click to Change Volume' will appear on the screen.
- Press the push button. The volume displayed on the screen will start blinking.
- Turn the push button clockwise to decrease volume or counterclockwise to increase volume
- Press the push button once when finished to lock the volume.
- Select a suitable tip for the pipette.
- PIPETMAN® M Connected is now ready to pipette.

Use the Menu Button

- Press the menu button to access the menu and adjust pipetting modes, speed, configuration, and to quit.
- Turn the push button to choose one of the modes. Press the menu button to confirm the selection.

NOTE

Pipette speed is set to six by default, but both the aspirate and dispense speeds can be adjusted in the speed menu.

Available Pipetting Modes

PIPET

Standard pipetting mode which aspirates and dispenses the volume set.

REPETITIVE

The pipette distributes the same volume repeatedly in an equal number of predefined aliquots.

MIX

Standard pipette mode followed by a mixing phase, i.e., aspirating and dispensing are repeated.

REVERSE

A preset amount of liquid is added during aspiration. After delivery, the excess volume remains in the tip and should be discarded.

CUSTOM

Enables up to 10 personalized pipetting protocol creation on a PC or an Android® tablet for transfer to the pipette. To use this mode, download the software from www.gilson.com/pipetmanmconnected or download the Android® app.

Turn Off the Pipette

Press and hold the push button for at least five seconds.

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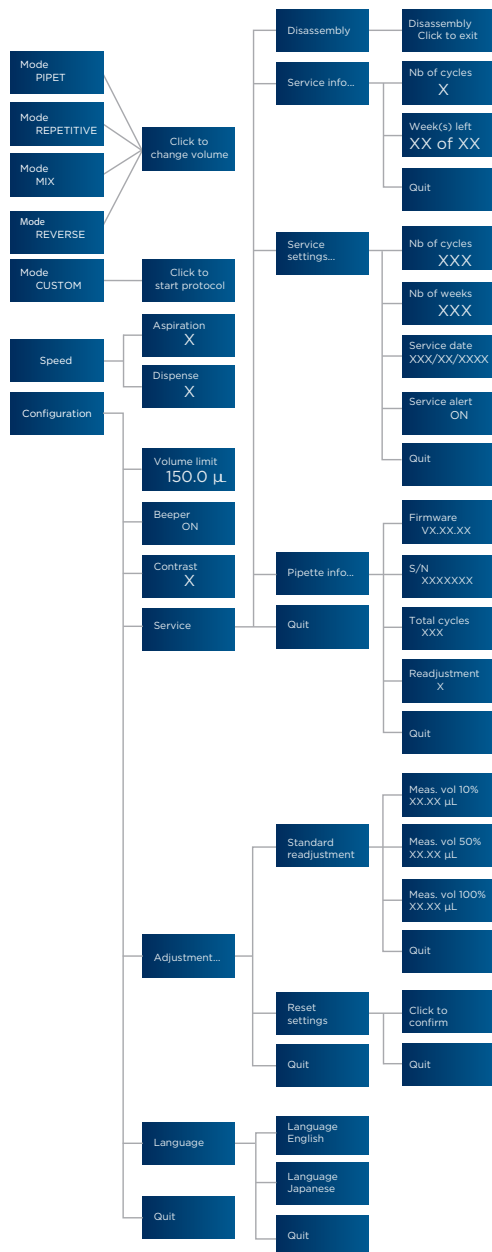
Menu Options

Bluetooth Connection

PIPETMAN M Connected can be connected to the PIPETMAN M Software and the Gilson apps either with the USB cable or using the Bluetooth® connection. To turn on Bluetooth, press the menu button, then press and hold the push button for 2-3 seconds until a blue light flashes.

Associated with TRACKMAN® Connected and its PipettePilot application, PIPETMAN M Connected will reduce risks of human errors by transferring verifiable pipetting parameters and commands in the pipette.

Please visit www.gilsonconnect.com to learn more about Gilson applications.

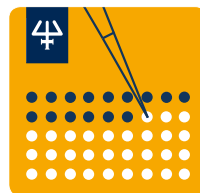
Figure 5
PipettePilot icon

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Gilson certifies on its sole responsibility that
PIPETMAN M Connected complies with the
requirements of the following European Directives:

2014/30/EU	Electromagnetic compatibility, EMC
2014/35/EU	Low Voltage Directive, LVD
2014/53/EU	Radio Equipment Directive

This Bluetooth enabled device also complies with the
following requirements:

USA, User information: Contains FCC ID: 2AAQS-ISP1507

Canada, User information: Contains IC: 11306A-ISP1507

Japan, TELEC certification n°207-16ISP5

The WEEE symbol (crossed-out wheeled
bin), according to the European Directive
2012/19/EU, indicates separate collection
for WEEE – Waste of Electrical and
Electronic Equipment.

Do not dispose electronic devices and
their batteries in a household bin, use the
recycling path in place in your country.

www.gilson.com/contactus

LT801595/A | ©2018 Gilson SAS. All rights reserved.